

Tamoxifen Protocol

- Tamoxifen
 - MP Biomedicals cat # 02156738
- Prepare fresh for injection
 - Measure out 10mg of Tamoxifen in 1.5ml eppendorf
 - Add 100 μ l 100% Ethanol
 - Vortex
 - Add 900 μ l Sunflower Seed Oil (Sigma-Aldrich cat # S5007)
 - Vortex
 - To ensure dissolved, heat 95°C for 1 min (waterbath)
 - Vortex to get a clear solution
- 100 μ l IP injection for 3 consecutive days