

Embryo Anesthetic

Flavio's recipe:

For 500ml:

5ml 10X MMR (final concentration is 0.1X)

0.1g 3-aminobenzoic acid ethyl ester (final concentration is 0.02%)

pH to 7.5 using a very small amount of NaOH.

Jen's Recipe:

For 100ml:

1ml 10 MMR (final concentration is 0.1X)

0.02g 3-aminobenzoic acid ethyl ester (final concentration is 0.02%)

0.02g NaBicarb

This should be around pH 7.0